

	Document #:	W-SW-0062	Rev:	A
	Title:	Simple Plex Explorer Version 5.0.1 Release Notes		
	Author:	Fred Westphal		
	Department:	Software		

1 SOFTWARE RELEASE NOTES



Software: Simple Plex Explorer

Version: 5.0.1

Released: July 2026

2 MINIMUM PC REQUIREMENTS

Operating System	Windows 11 64-bit
Processor	Core i7 – 6 th Generation
Memory	16 GB
Storage	500 GB
Display Resolution	1600x900 (with 100% Display Scaling), 1920x1080 (with 125% Display Scaling)
USB	1 Ports – USB 2.0 Type A (1 for barcode scanner)

3 INSTALLATION

1. Uninstall previous version of Simple Plex Explorer (if present)
 - a. Open Settings – Apps – Installed apps
 - b. Click on the ‘...’ for the existing Simple Plex Explorer app, select Uninstall and follow the prompts.
2. Install new version of Simple Plex Explorer
 - a. Double click on the installation file, Simple Plex Explorer 5.0.1.x.msi and follow the prompts.

4 COMPATIBILITY WITH OTHER SOFTWARE

N/A

5 VERSION 5.0.1.6 SIMPLE PLEX EXPLORER

5.1 FEATURES

N/A

	Document #:	W-SW-0062	Rev:	A
	Title:	Simple Plex Explorer Version 5.0.1 Release Notes		
	Author:	Fred Westphal		
	Department:	Software		

5.2 IMPROVEMENTS AND FIXES

1. FIX(REX-327): For 21 CFR Part 11 mode, resolves login issues in some domain-based Active Directory scenarios where organizational policy disallows password-based Windows sign-in (for example, when Windows Hello is enforced). A “Login failed” message was shown despite a valid username and password being entered.
2. FIX(REX-333): For 'Single Analyte QC Report 1' export, resolves issue where the 'Ella Analyzer Description' field in the first section was incorrectly showing analyzer name.
3. CHANGE(REX-331): Minor changes to PDF exports (table splitting behavior to reduce gaps, etc.)

5.3 KNOWN ISSUES

N/A

6 REVISION HISTORY

Table 6-1. Revision history

Revision	ECO	Change Descriptions	Initiated by
A	ECO-W-006182	Initial revision	Fred Westphal