

Compass for SW 6.1.0 and Embedded Software Installation Guide: Steps to Update Your Jess System

Jan 18, 2022

Note: *Both the Jess instrument (embedded) software and Compass for Simple Western **must be updated to enable** all the new **Stellar Assay** features. It is recommended that you update Compass software prior to updating the embedded software.*

Downloading Software (With an Internet Connection)

Downloading Compass for SW 6.1.0

If the instrument PC is connected to the internet, open Compass for Simple Western select **Help > Check for Updates** to download the latest version of the software. Follow the on-screen instructions to complete the software installation.

Downloading Embedded Software

If the instrument PC is connected to the internet, then the update can be completed by having Compass for Simple Western update the instrument software from the latest version on our website. To check for embedded updates, you will first need to make sure that the Compass for Simple Western software is connected to the instrument by selecting the **Instrument** menu, then **Connect**. Once connected, select **Instrument** in the main menu, then select **Update** and **Network**.

Updating the Software with No Internet Connection

Updating Compass for Simple Western 6.1.0

If the instrument PC is not connected to the internet, follow these steps:

- 1) Visit <http://www.proteinsimple.com/compass/downloads/> to download the installer using a PC with an internet connection.
- 2) Transfer the Compass software installer to a USB drive.
- 3) Put the USB drive in a USB slot on the Jess computer.
- 4) Run the installer.

Updating Embedded Software (No Internet Connection)

If the instrument PC is not connected to the internet, follow these steps:

- 1) Contact Technical Support or your local Field Applications Specialist to receive the latest embedded installer file.
- 2) Transfer the embedded installer file (*embedded-rg-2.1.2021.09.01.00.26.14.bc9a11c63.tar.gz*) to a USB drive.
- 3) Put the USB drive in a USB slot on the Jess computer.
- 4) Make sure that the Compass for Simple Western software is connected to the instrument by selecting the **Instrument** menu, then selecting **Connect**.
- 5) Select the **Instrument** menu, select **Update**, then select **Local...**
- 6) In the open dialog window, browse to the location of the embedded installer file (*embedded-rg-2.1.2021.09.01.00.26.14.bc9a11c63.tar.gz*), select the file and click **Open** to initiate the update of embedded software.