



Release Notes – Compass for SW software

Version 7.1.0

P/N 805-0003, Revision G

Released November 2025

Modified release notes from Rev F to enhance clarity regarding existing functionality; Rev F and G include same software changes across Leo, Jess and Abby software applications.

Recommended PC requirements for Leo

Operating System	Windows 11 Pro/Enterprise 64 bit
Processor	Intel Core i7 or better
Memory	16 GB
Free Disk Space	100GB
Ethernet Ports	2 (1st port for Instrument, 2nd port for network access)

Recommended PC requirements for Jess/Abby

Operating System	Windows 11 Pro/Enterprise 64 bit
Processor	Intel Core i3 or better
Memory	8 GB or more
Free Disk Space	100GB
Ethernet Ports	2 (1st port for Instrument, 2nd port for network access)

Installation

Open the installer file to start the installation.

NOTE: User can download the installer from Bio-Techne download center website via 'Help | Check for Updates' feature in Compass software or by visiting the link: [Bio-Techne | Compass for Simple Western Download Center](#). Please note that the direct software update from version 3.x to version 4.0.0 and higher is no longer supported. User will need to download installer and perform a full installation.

NOTE: If you are using Peggy Sue, Sally Sue or NP1000, Compass for Simple Western version 6.0 and higher does not support these instruments. You can still use software versions 5.x and lower to create, load and run assays on these instruments, and to analyze your data.

NOTE: If you are using Wes, Compass for Simple Western version 7.0 and higher does not support the instruments. You can still use software versions 6.3 and lower to create, load and run assays on these instruments, and to analyze your data.

Compatibility with Instrument software and Authorization Server

Following software is tested and released with Compass for SW 7.1.0.

Instrument Software:

Leo Instrument software 1.2, Jess 2.2 and Abby 1.4 to support respective instruments.

Compass Authorization Server for SW: 3.0.0

Compass for SW 6.2.0 or above users are required to upgrade to Compass Authorization Server for SW version 3.0.0 (if not already) to establish secure connection between Compass for SW and Authentication server (CAS).

If Compass for SW tries to connect to older version of CAS (below version 3.0.0), a warning is displayed that Compass Authorization Server is out of date and is connecting in insecure manner.

For more details on Compass Authorization Server, contact Technical Support.

7.1.0 Compass for SW

Features

1. Added Fluorescence support for Leo instrument (CSW-1656).
Note: For Fluorescence assay runs, instruments with Leo 1.0 require hardware upgrade and instrument software Leo 1.2. If the upgrade is not done, users can still perform Chemiluminescence runs with Compass for SW 7.1 and Leo 1.0.
2. Compass for SW 7.1 restricts users to upgrade 'Leo instrument software from 1.0 version without hardware upgrades (CSW-2038)
Note: Contact ProteinSimple Technical Support for assistance on how to obtain the latest update. Jess and Abby embedded upgrades from network shall work fine.
3. Added support for new sample plate for Leo where row N is trough without well (CSW-1919)
Note: Compass for SW 7.1 will display the new sample plate image in Layout view, even for run files generated with Compass for SW 7.0
4. Compass for SW shall allow users to freely set Primary antibody incubation time instead of pre-selecting values from a drop down(CSW-1984)
5. Provided an option 'Well Info from' to choose Sample or Primary or Secondary Antibody name to load default concentration values in Standard Curve analysis settings (CSW-1841)
Note: The rule of concentration applied will remain the same as in v7.0
6. Removed Jess-Stellar assay support from Compass for SW (CSW-1619)

Improvements and Fixes

General

1. Assay Template changes made during the run are now saved after the run completes (19050 or CSW-306)
2. Increased Standard trace line thickness to differentiate between Sample and Standard when overlapped in graph view (CSW-1914)

3. Fixed the issue where copy-paste from excel to Compass Assay template caused an error and resulted in only a single row being pasted (CSW-1986)
4. Corrected 'Corr. Area' calculation when the Peak Group does not contain reference caps from all 4 cartridges and selected Corr. Cart is Cartridge Average (CSW-1988)
5. Corrected order of groups in Peaks table as per the cartridge number when groups are created per cartridge (CSW-1997)
6. Improved lane contrast for independent cartridge and probe scale selection with or without Baseline On (CSW-2026)
7. Fixed an issue with Jess Chemi PN assay where ladder channel was set to NIR channel by default. Now it is set to CHEMI as default (CSW-2031)
8. Fixed an issue in the Leo Assay Setup Wizard where the number of sample rows would reset to match the number of cartridges selected (CSW-1929)
9. Fixed an issue in Lane View, Capillaries and Capillaries Groups table where caps for non-ladder cartridges were missing when the ladder was selected from a specific cartridge instead of being assigned per cartridge in analysis ladder settings (CSW-2096, CSW-2104, CSW-2105)
10. Fixed an issue in the selected Graph View where the 'Name Peak' option was missing from the right-click context menu for the caps of non-ladder cartridges when the ladder was selected from a specific cartridge in the Analysis Ladder settings (CSW-2103)
11. Fixed an issue in export tables where Peaks.txt file has missing peaks for the caps of non-ladder cartridges when the ladder was selected from a specific cartridge in the Analysis Ladder settings (CSW-2107)
12. Fixed an issue in the History log where changes to "Independent Probe Scale" in lane options were not being recorded (CSW-2123)
13. Fixed an issue where the Peak Area for Probe 2 was not updating in the Peak table and Graph view when the Area Calculation Analysis setting was set to 'Dropped Lines (manual)' (CSW-2140)
14. Fixed an issue where Lane contrast adjustments were not preserved after reopening a saved run file (CSW-2141)
15. Jess flat field specs are updated for manufacturing to account for expected component batch variations (CSW-2037)

Report

1. Added the channel name to the normalization analysis settings for Leo and Jess in the run reports (CSW-2007)

Known Issues

General

1. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
2. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
3. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)

4. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
5. Start Assay and Blot dialogs can be displayed at the same time. (17467)
6. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
7. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
8. Channel Column appears for Peak Groups in Abby run file Reports (CSW-50 or 21341)
9. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (CSW-1718 or 21394)
10. Lane View – Probe 2 Ladder Lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (CSW-55 or 21410)
11. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
12. When Exposure settings contain decimal values, folder name created from 'Export Tables' action in Compass does not display decimals (21613)
13. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)
14. Corrected area column is missing when Normalization is enabled and when ladder cap is set as reference cap (21676 or CSW-567)
15. For Locale Chinese and Japanese, in assay perspective well name changes are not retained after Compass restarts. Report and export tables show empty well names (21677)
16. In graphs view, when 'stack plots' view is on, consecutive zooming in multi-lane view resets initial zoom action on the Y-axis to display default scale. (CSW-289)
17. Compass crashes with an internal error dialog when zoom action is performed repeatedly in the graph view in an empty graph area (CSW-313)
Solution: Close the error dialog and launch Compass for SW again, there is no impact on run result data or analysis settings due to this issue.
18. Export of analysis settings prompts to save in a path different from the run file path (CSW-317)
Solution: One can navigate to intended path and save the analysis settings
19. Independent Probe Scale option is checked in Lane View even after the setting was unchecked in Graph View (CSW-321)
20. In images view, capillary selection on Probe 2 images disappears when navigated between standards and samples views (CSW-325)
21. History log incorrectly displays standard curve name without channel information although peak name assigned to standard curve is not changed (CSW-583)
22. Assay perspective, Edit->Cut option is always disabled even when text is selected (CSW-1715)
Solution: Ctrl + X operation works on text
23. Group plots are not available for Leo runs (CSW-929).
Note: For Jess and Abby runs, no change in Group Plots calculation. There are plotted for Sample Names and Primary Names only and are not changed per Name and Attribute selections in Analysis settings > Groups selection.

24. In rare cases, when a position of standard is changed and saved in run file, it may not show the same position when run file is opened again (CSW-1725)
25. If a channel has no peaks, the run report can potentially include blank pages (CSW-2077)
26. Report does not display the standard curves with error (CSW-2083)
27. While the run is in progress, the ladder channel may not populate in the Analysis settings (CSW-2118)
Note: The ladder channel populates as expected after the run completes
28. Compass displays an error dialog when "Set to PC Time" is executed while the instrument is in an error state (CSW-2112)
Solution: Clear/reset the reported error in the instrument before performing "Set to PC Time"
29. Undo and Redo actions do not reflect changes when the Analysis settings window is open (CSW-2111)
Solution: To ensure Undo/Redo functionality works correctly, close the Analysis settings window before performing Undo or Redo actions
30. Copying data from Excel to the Compass assay template will omit the last row if that row has not been edited in Excel. As a result, the last row might not appear in the Compass assay template view after pasting. This usually happens when initial excel contents were copied from CSV file (CSW-2157)
Solution: Edit the last row in Excel before copying, or manually add the missing row in Compass after pasting.
31. Standard curves and cartridge corrections will fail to load correctly if peak names include channel identifiers that software application uses, such as '(CHEMI)', '(NIR)', or '(IR)'—for example, 'Peak 1 (CHEMI)' or 'Peak 2 (NIR)' (CSW-2163)
32. The pdf run report Peaks table shows truncated data if it contains more than 18 columns (CSW-2166)
33. If a peak group has multiple peaks with the same name but different channels, and those peaks are used with standard curves and cartridge corrections, Compass will not load the standard curves or cartridge corrections properly (CSW-2171)
Solution: Provide unique peak names or create separate peak groups and peak names to use with standard curves and cartridge corrections for Compass to load standard curves or cartridge corrections properly.
34. In few cases of log-log standard curve fit in Compass for SW, R-squared value returns negative or incorrect reading due to an issue in the R^2 formula.
Solution: Data point display and Log-Log curve fit of your standard curve is correct except for R-squared formula, you can export data and manually calculate R-squared value.

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

1. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)

2. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272 or CSW-1072)
3. Cannot copy text to Windows clipboard from Notes or Labels (21289 or CSW-1657)
4. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
5. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
6. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
7. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
8. Underlines are not aligned for angled and wrapped text Lane labels (21353)
9. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)
10. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)
Solution: Open the exported .tiff image with 'Photos' application in Windows and then use right-click option to copy the image into Word/PowerPoint/etc.
11. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)
12. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly
Solution: Perform Re-crop to update the Figure Band Label displays (21440)
13. Compass freezes when trying to generate a BMP format for the Replex TP Lane view image when both probe 1 and 2 are selected (CSW-2082)

RePlex™

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
3. Selecting Independent Probe Scales for both Probe 1 & Probe 2 results in vertical lines for Probe 2 in Graph view (CSW-2133)
Solution: Enable Auto-scale and disable it again to fix this issue.

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability
Solution: Use the same fixed contrast levels for added runs (16226)
3. Lane order resets when user switches View from Single to All (21270)
4. Lane order resets after user changes Lane Options (21271)

History

1. Lane View Options and Graph view option changes are not saved correctly in History log (21443)
Solution: Pre-existing access-controlled Jess non-Replex assays that were created with versions earlier than 6.1 will have this issue. If history log entries are critical to the system, perform an initial save of the assay file. Logs of any changes after the first save will be displayed correctly.

1.2 Leo Instrument Software

Features

1. Release of instrument software for Leo to enable Fluorescence features (ESW-265).
2. Added additional instrument self-tests to verify functionality for fluorescence related features (ESW-1344, ESW-1345)
Note: Contact ProteinSimple Technical Support for assistance on how to upgrade from 1.0 to 1.2 instrument software.

Improvements and Fixes

1. Modified self-test algorithms and specifications to improve reliability and optimize self-test duration (ESW-1329, ESW-1340, ESW-1423, ESW-1424).
2. Fixed inconsistencies with the movie camera rotation self-test results (ESW-1444).

Known Issues

N/A

2.2 Jess Instrument Software

Features

1. None

Improvements and Fixes

1. Fixed an issue in the low vacuum calibration test reporting, which previously under-reported the difference between measured values by a factor of four. The calculation now accurately reflects the true difference, ensuring correct pass/fail reporting in Jess self-test reports (ESW-97)
2. Updated the specification for the FL LED difference in on/off states during self-test. The new specification aligns with the intended distinction between manufacturing, field service, and customer usage, allowing for more accurate assessment of LED degradation over time (ESW-1260)

3. Extended the expiration window for the HPLED Calibration tool from 6 months to 12 months, reducing the frequency of required recalibration and improving maintenance efficiency (ESW-1357)
4. Updated the vacuum leak self-test user specifications to 0.2 in Hg/min and service specification to 0.1 in Hg/min. This is to remove a false failure of vacuum leak self -test while system is still performing to the intended specifications (ESW-1396)
5. Fixed this: When a run is performed with chemi detection setting other than HDR in the detection profile of Jess assay, chemi channel will not show any data (ESW-142).
Solution: Select High Dynamic Range (HDR) as the Chemi Detection profile to perform chemi based runs on Jess. Details of HDR can be found in Compass for SW user guide
6. Fixed an issue where the filter wheel could remain on a multi-fluor position after acquiring chemi images. The filter wheel is now explicitly moved to a position for chemi imaging, ensuring proper operation even after previous multi-fluor acquisitions (ESW-142)

Known Issues

1. On rare occasions when attempting to close the door, the instrument will not properly recognize the hand sensor input and close (17590)

1.4 Abby Instrument Software

NOTE: Requires Compass for SW version 6.0 or higher

Features

1. None

Improvements and Fixes

1. Fixed an issue in the low vacuum calibration test reporting, which previously under-reported the difference between measured values by a factor of four. The calculation now accurately reflects the true difference, ensuring correct pass/fail reporting in self-test reports (ESW-1288)
2. Updated the camera aperture self-test specification from $\pm 15\%$ to $-15\%/+20\%$. This change provides a more accurate and appropriate tolerance for the Abby platform, with no changes made to Jess (ESW-1261)
3. Fixed a bug in the Lamp Set Point (calibration) self-test where the reported percentage difference was incorrect, potentially allowing a failing result to be marked as a pass. The calculation now correctly reflects the actual change, ensuring accurate test results (ESW-1388)
4. Updated the vacuum leak self-test user specifications to 0.2 in Hg/min and service specification to 0.1 in Hg/min. This is to remove a false failure of vacuum leak self -test while system is still performing to the intended specifications (ESW-1396)

Known Issues

1. None

7.0.0 Compass for SW

March 2025

Features

1. Added support for Leo instrument.
2. Added ability to perform group edits in assay template view and made more usability improvements. Compass for SW now supports split well content for names and attributes, copy-paste operations and key navigation with up/down keys. It also allows opening edit text box with Enter key etc., in the template view (CSW-411)
3. In the Added runs scenarios, added sub menus to perform file actions such as Close, Save As, Export Tables, Export Spectra – Text and ANDI Formats and Run Reports on user selected run file (CSW-960, CSW-961, CSW-1243)
4. In the Added runs scenario, under the File-> Export Tables menu, added 'Batch Exports' submenu option to let users create a table that combines the data of all exported files (CSW-1278)
5. Added a toggle button to enable or disable Grouping Across Runs for Added runs scenarios (CSW-955)
6. Added ability to choose between linear or log axes scales for standard curves for both x and y axes (CSW-349)
7. Added a new fit type "Log-Log" for Standard Curve (CSW-1091)
8. Added Weighting factor support for all fit types in Standard Curve (CSW-1092)
9. Added a new "Std. Curve" tab/table to show curve fit results in the analysis view (CSW-966). Additional parameters such as Min (a), Max (d), EC50 (c), Y0 (c), Hill Slope (b) and Slope (m) are displayed now in Standard Curve table (CSW-1313)
10. In standard curve graph view, added legend to show std curve formula, R-squared value and other key parameters, standards are displayed as blue dots and samples in red (CSW-1114)
11. Standard Curve analysis settings now display cap locations, Sample names and attributes to help user to select standard caps and set known concentrations accordingly (CSW-773)
12. Introduced new column "Known Conc." to display the concentrations entered in Analysis settings for Standard Curve (CSW-740)
Note: The existing column "Conc." will display the calculated concentration for all caps in the peak group along with the Standard capillaries. Hence, the concentration data for Standard capillaries will be different from the data in 6.3 or lower Compass SW versions.
13. Added a new page under Analysis Settings for Grouping. Grouping can now be done by sample, primary or secondary names or attributes. In previous versions of Compass Grouping option was under Edit->Preferences menu (CSW-406, CSW-825)
Note: Compass for SW 7.0 will display different data for Grouping by default for Jess runs as 'Secondary Name' is not selected in Analysis Group Settings. However, user can change Grouping preferences and can match the way it was in 6.3 and lower versions.
14. Introduced new help menu "Online Antibody Database" to access Simple Western Antibody Database (CSW-716)

15. Removed Twitter feature from Compass for SW under Preferences (CSW-626)
16. Warnings and errors from Leo instrument are now displayed in the Instrument status or run progress bar area of Compass for quicker issue visibility (CSW-1673)

Improvements and Fixes

General

1. Fixed the issue of assay Template changes. Changes made during the run are now saved after the run completes (19050 or CSW-306)
2. Added import and export buttons in the assay template view for better visibility. These operations can also be performed from the File menu (CSW-480)
3. Improved visibility of e-grams in graph view for better view in semi bright settings by making the lines thicker (CSW-312)
4. Fixed an issue where 'zoom out' only worked properly for first standard curve graph. Also, when reanalysis of standard curve happens in a custom layout, a new standard curve was generated while old one is still displayed, fixed this scenario (21039 or CSW-704, CSW-1083)
5. Improved history logging. In previous versions of Compass for SW, detailed log entries were computed by analyzing differences in edit.xml files. This method is difficult to maintain as we continue to add more functionality. Therefore, changed the logging mechanism to track setting changes by user action and generate detailed log entries from those changes as they happen. Detailed log entries for the most part will appear same as in previous versions of Compass for SW. No information is lost but changes are consolidated (CSW-1654)
6. Data difference: In compass for SW 6.2 or above, for Replex runs, when Force manual standard operation is performed on one probe, it did not reflect for another, fixed this issue (CSW-1547)
7. Fixed the issue where hidden cap was displayed in Lane view when it's hidden from experiment tab view (CSW-1527)
8. Fixed an NPE error that was causing a non-Replex assay to not open or be visible to user when a Replex assay was open in Compass (CSW-694)
9. For Jess and Abby, moved 'Group by Probe' option from Preferences to Analysis Settings and user selection of the view is saved in the run file (CSW-494)
10. In Jess, fixed the issue of inaction when 'Remove All' button is selected in Chemi detection profile window (CSW-625)
11. Signal to Noise settings – validation checks are added to ensure Start < End (21229 or CSW-270)
12. When Custom noise region – start and end values are changed in Signal to Noise setting is changed, in Compass for SW 6.2 or above, the end region was not displayed correctly in the graph view, fixed this issue (CSW-1056)
13. Compass in access control mode cannot connect to an instrument running an assay in non-access control mode and vice-versa (CSW-1545)
14. Annotations tab was displayed in analysis view in CFR mode, fixed it as annotations are not supported in CFR mode (CSW-1640)
15. Changed the default colors of NIR/IR Channels in probe 2 of Replex runs to improve visualization in Graph view (CSW- 941)
16. Fixed issue with lane selection for RePlex – Probe 2 lane is now selected if channel added after capillary selection done in Lane View (21388 or CSW-1625)

17. Issue with standards curve graph tab not displayed after undo redo actions in analysis settings is fixed (CSW-72)
18. Fixed the issue with custom contrast changes being ignored if independent probe scale was not selected. Custom contrast changes are now being reflected in the annotation regardless of the independent probe scale (21455 or CSW-1400)
19. Fixed issue of peak name group entries not logged (21692 or CSW-1629)
20. Fixed an issue to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected (CSW-1626 or 19434)
21. Groups with same peak names are now displayed with respective channel info. If there are peaks with same name in different channels for Compass SW v6.3, then previous channel data was getting overwritten with last analyzed channel (CSW-977)
22. Fixed an issue with baseline subtraction with Interpolate baseline type and Area calculation set to Dropped Lines or Dropped lines manual. This change can show different peak height and width data between Compass version 6.3 and 7.0. However, peak area remains same (CSW-1225)
23. Fixed Area Corrections for caps with multiple loading controls to consider the reference caps from multiple loading controls and not just the last loading control (CSW-1222)
24. Run files with normalization enabled uses peak corrected area for standard curve, Y-axis label is corrected to show peak corrected area in such cases (CSW-1809)
25. Export spectra in text format now contain data for standard curves if they are present in run file (CSW-912)
26. Fixed the issue of Compass for SW where non-digital signature actions are not being logged in the CAS audit trail due to session expiration between Compass and Compass Authentication Server (CSW-1310).
Note: The session expiration time between Compass and CAS is independent of the Compass screen auto-lock inactivity time that the user can set on Compass user interface.
27. Fixed the CAS disable dialog box workflow issue. When Compass is launched in Access Control mode and it fails to connect with CAS, the error dialog will appear without the 'Disable' option unless Compass was started with 'Run as administrator' (CSW-940)

Known Issues

General

1. Add Run - Scale incorrect when run files of different sizes are added. (18915)
2. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
3. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
4. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
5. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
6. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)

7. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
8. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
9. Start Assay and Blot dialogs can be displayed at the same time. (17467)
10. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
11. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
12. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
13. Standard Curve tab disappears after user Adds Run without Std Curve and reanalysis is triggered on second run (21322)
14. Channel Column appears for Peak Groups in Abby run file Reports (21341)
15. Standards Graphs - Exposures in legends should be in black when the plots are stacked (21351)
16. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (21394)
17. Lane View – Probe 2 Ladder Lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (21410)
18. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
19. When Exposure settings contain decimal values, folder name created from 'Export Tables' action in Compass does not display decimals (21613).
20. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)
21. Corrected area column is missing when Normalization is enabled and when ladder cap is set as reference cap (21676 or CSW-567)
22. Add runs – 'save', 'close' and 'close all' operations
 - Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed file. The user can then save and close.
 - Selecting File->Close All does not close all the files after saving (19048)
 - When two runs are open, selecting File->Close on a specific run will close the other run (21797).
23. For Locale Chinese and Japanese, in assay perspective well name changes are not retained after Compass restarts. Report and export tables show empty well names (21677)
24. In graphs view, when 'stack plots' view is on, consecutive zooming in multi-lane view resets initial zoom action on the Y-axis to display default scale. (CSW-289)
25. Compass crashes with an internal error dialog when zoom action is performed repeatedly in the graph view in an empty graph area (CSW-313)
Solution: Close the error dialog and launch Compass for SW again, there is no impact on run result data or analysis settings due to this issue.
26. Export of analysis settings prompts to save in a path different from the run file path (CSW-317)

Solution: One can navigate to intended path and save the analysis settings

27. Independent Probe Scale option is checked in Lane View even after the setting was unchecked in Graph View (CSW-321)
28. In images view, capillary selection on Probe 2 images disappears when navigated between standards and samples views (CSW-325)
29. History log incorrectly displays standard curve name without channel information although peak name assigned to standard curve is not changed (CSW-583)
30. Assay perspective, Edit->Cut option is always disabled even when text is selected (CSW-1715)

Solution: Ctrl + X operation works on text

31. Group plots are not available for Leo runs (CSW-929).

Note: For Jess and Abby runs, no change in Group Plots calculation. There are plotted for Sample Names and Primary Names only and are not changed per Name and Attribute selections in Analysis settings > Groups selection.

32. In rare cases, when a position of standard is changed and saved in run file, it may not show the same position when run file is opened again (CSW-1725)

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

1. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)
2. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272)
3. Cannot copy text to Windows clipboard from Notes or Labels (21289 or CSW-1657)
4. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
5. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
6. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
7. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
8. Underlines are not aligned for angled and wrapped text Lane labels (21353)
9. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)
10. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)
11. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)
12. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly

Solution: Perform Re-crop to update the Figure Band Label displays (21440)

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability

Solution: Use the same fixed contrast levels for added runs (16226)

3. Lane order resets when user switches View from Single to All (21270)
4. Lane order resets after user changes Lane Options (21271)

History

1. Lane View Options and Graph view option changes are not saved correctly in History log (21443)
Solution: Pre-existing access-controlled Jess non-Replex assays that were created with versions earlier than 6.1 will have this issue. If history log entries are critical to the system, perform an initial save of the assay file. Logs of any changes after the first save will be displayed correctly.

1.0 Leo Instrument Software

Features

Initial release of instrument software for Leo

Improvements and Fixes

N/A

Known Issues

N/A

6.3.0 Compass for SW

August 2023

Features

1. Added the ability to manually adjust the start and end pixel position in analysis settings for Protein Normalization baseline (CSW-417).

Improvements and Fixes

General

1. Compass for SW and Instrument user guides offline copies are no longer installed by default with Compass for SW installer. Latest user guides can be obtained from Bio-Techne website (CSW-425)
2. Added informational message in instrument properties window to restart compass and instrument after selecting 'Set to PC time' button (CSW-426)
3. Added a reminder in the Start UI dialog to fill all empty wells in rows that are being used in the assay before starting a run (CSW-520)
4. Fixed issue with incorrect incubation wash steps generated by Compass for SW for Stellar assays. This issue occurs when new assay is created from File->New Assay Menu and the fix will not impact performance of runs already performed (CSW-457)
5. Fixed the issue of the inability to select and copy multiple caps in Capillaries tab under 'View Selected' option (CSW-472, CSW-482)
6. Fixed the annotation issue with probe 1 channel (s) selected for RePlex assays. Annotation can be done with either probe selection, 'index out of bounds' error no longer appears (CSW-353)
7. Fixed the issue where IV plots are not exported into run report for certain runs (CSW-354)
8. Fixed issue with contrast intensity of the ladder cap in lane view. Ladder cap intensity is now independent of rest of the caps and is set per peak fit view range in analysis settings, sample lane contrast is set per peak fit view range in analysis settings as well (CSW-499, CSW-500)
9. Normalization settings of Stellar Total Protein runs are now added in the run report (CSW-498)
10. Fixed the issue of Compass for SW displaying 'Cannot connect to server' warning dialog even when authentication server was connected (CSW-538).

Known Issues

General

1. IV Plot title for Jess run displays "Cycle 1", instead of being blank (18962)
2. History Log - Manual standard changes log message is not easily understandable. (18882)
3. Add Run - Scale incorrect when run files of different sizes are added. (18915)
4. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
5. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
6. Analysis Settings for Peak Names -> Standard Curves -> Concentration - units of precision do not match report. (18905)

7. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
8. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
9. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)
10. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
11. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
12. Start Assay and Blot dialogs can be displayed at the same time. (17467)
13. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
14. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
15. Standard Curve tab disappears upon reanalysis of data if user selects to Cancel settings dialog
Solution: Go to Edit > Analysis... > Standard Curve > OK (19012)
16. Assay Template changes made during the run are not saved after the run completes (19050 or CSW-306)
17. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
18. Zoom out only works properly for first standard curve graph (21039)
19. Signal to Noise settings – validation checks do not exist to ensure Start < End (21229)
20. File > Close All does not close all run files if one of them needs to be saved (21320)
21. Standard Curve tab disappears after user Adds Run without Std Curve and reanalysis is triggered on second run (21322)
22. Channel Column appears for Peak Groups in Abby run file Reports (21341)
23. Standards Graphs - Exposures in legends should be in black when the plots are stacked (21351)
24. Lane selection for RePlex – Probe 2 lanes not selected if channel added after capillary selection done in Lane View (21388)
25. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (21394)
26. Lane View – Probe 2 Ladder Lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (21410)
27. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
28. When Exposure settings contain decimal values, folder name created from 'Export Tables' action in Compass does not display decimals (21613).
29. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)
30. Corrected area column is missing when Normalization is enabled and when ladder cap is set as reference cap (21676 or CSW-567)
31. Add runs – 'save', 'close' and 'close all' operations

- Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed file. The user can then save and close.
 - Selecting File->Close All does not close all the files after saving (19048)
 - When two runs are open, selecting File->Close on a specific run will close the other run (21797).
32. For Locale Chinese and Japanese, in assay perspective well name changes are not retained after Compass restarts. Report and export tables show empty well names (21677)
 33. Standards curve graph tab does not appear after an undo-redo action (CSW-72)
Solution: Standards curve tab can be restored by opening analysis setting and clicking on ok button.
 34. In graphs view, when 'stack plots' view is on, consecutive zooming in multi-lane view resets initial zoom action on the Y-axis to display default scale. (CSW-289)
 35. Compass crashes with an internal error dialog when zoom action is performed repeatedly in the graph view in an empty graph area (CSW-313)
Solution: Close the error dialog and launch Compass for SW again, there is no impact on run result data or analysis settings due to this issue.
 36. Export of analysis settings prompts to save in a path different from the run file path (CSW-317)
Solution: One can navigate to intended path and save the analysis settings
 37. Independent Probe Scale option is checked in Lane View even after the setting was unchecked in Graph View (CSW-321)
 38. In images view, capillary selection on Probe 2 images disappears when navigated between standards and samples views (CSW-325)
 39. History log incorrectly displays standard curve name without channel information although peak name assigned to standard curve is not changed (CSW-583)

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

13. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)
14. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272)
15. Cannot copy text to Windows clipboard from Notes or Labels (21289)
16. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
17. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
18. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
19. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
20. Underlines are not aligned for angled and wrapped text Lane labels (21353)
21. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)

22. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)
23. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)
24. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly
Solution: Perform Re-crop to update the Figure Band Label displays (21440)
25. For RePlex runs when "Independent Probe Scale" setting is Off and the user modifies the Contrast in the Lane View, the newly created Figures don't have the same contrast
Solution: Use Custom Contrast in Crop Regions to adjust contrast as desired (21455)

RePlex™

3. Plot colors for RePlex runs may not be in the order expected. (18943)
4. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
5. Unable to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected.
Solution: Temporarily unselect the Total Protein channel, then the context menu appears, allowing the user to make any necessary changes (19434)

Lane

5. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
6. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability
Solution: Use the same fixed contrast levels for added runs (16226)
7. Lane order resets when user switches View from Single to All (21270)
8. Lane order resets after user changes Lane Options (21271)

History

2. Lane View Options and Graph view option changes are not saved correctly in History log (21443)
Solution: Pre-existing access-controlled Jess non-Replex assays that were created with versions earlier than 6.1 will have this issue. If history log entries are critical to the system, perform an initial save of the assay file. Logs of any changes after the first save will be displayed correctly.
3. Removal of Peak name groups entries are not logged (21692).

6.2.0 Compass for SW

Released October 2022

Features

1. Added Manual Peak Integration feature. Peak area can now be adjusted by dragging start and end of the peak when 'Dropped Lines (manual)' option is selected under Peak Find analysis settings (CSW-1)
2. Added new 'Interpolate' option for Baseline type in Peak Fit analysis settings. When selected, Compass for SW lets you interpolate baselines from the base of each peak when chosen fit types are Dropped Lines or Dropped Lines (manual) (CSW-10)
3. Added ability to perform undo - redo actions within Compass for SW for assay and analysis views (CSW-92)
4. Added Ability to adjust peak table columns (peaks, capillaries, and experiment tables) in the analysis view (CSW-132).

Note: These column selections are limited to Compass views only, not reflected in exports and reports.

Improvements and Fixes

General

1. Help resources are moved to Bio-Techne website as proteinsimple.com site retired (CSW-80)
2. Removed unused Log4j library from code to avoid any vulnerability threats (CSW-93)
3. Bio-Techne Academy link is added in the help menu to provide easy access to learning portal (CSW-96)

RePlex™

1. In some Re-Plex standards run, Y-axis displays chemi units instead of Fluorescence (21429 or CSW-52)
2. Corrected default Peak Fit settings for Low MW assays so that they are the same for Abby and Jess, with or without Total Protein (CSW-74)
3. Fixed issue with missing Total Protein data after generating std curves in Jess RePlex TP run (CSW-450)

Known Issues

General

1. IV Plot title for Jess run displays "Cycle 1", instead of being blank (18962)
2. History Log - Manual standard changes log message is not easily understandable. (18882)
3. Add Run - Scale incorrect when run files of different sizes are added. (18915)
4. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
5. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
6. Analysis Settings for Peak Names -> Standard Curves -> Concentration - units of precision do not match report. (18905)
7. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
8. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)

9. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)
10. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
11. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
12. Start Assay and Blot dialogs can be displayed at the same time. (17467)
13. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
14. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
15. Standard Curve tab disappears upon reanalysis of data if user selects to Cancel settings dialog
Solution: Go to Edit > Analysis... > Standard Curve > OK (19012)
16. Assay Template changes made during the run are not saved after the run completes (19050 or CSW-306)
17. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
18. Zoom out only works properly for first standard curve graph (21039)
19. Signal to Noise settings – validation checks do not exist to ensure Start < End (21229)
20. File > Close All does not close all run files if one of them needs to be saved (21320)
21. Standard Curve tab disappears after user Adds Run without Std Curve and reanalysis is triggered on second run (21322)
22. Channel Column appears for Peak Groups in Abby run file Reports (21341)
23. Standards Graphs - Exposures in legends should be in black when the plots are stacked (21351)
24. Lane selection for RePlex – Probe 2 lanes not selected if channel added after capillary selection done in Lane View (21388)
25. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (21394)
26. Lane View – Probe 2 Ladder Lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (21410)
27. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
28. When Exposure settings contain decimal values, folder name created from 'Export Tables' action in Compass does not display decimals (21613).
29. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)
30. Corrected area column is missing when Normalization is enabled and when ladder cap is set as reference cap (21676)
31. Add runs – 'save', 'close' and 'close all' operations
 - Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed file. The user can then save and close.
 - Selecting File->Close All does not close all the files after saving (19048)

- When two runs are open, selecting File->Close on a specific run will close the other run (21797).
32. For Locale Chinese and Japanese, in assay perspective well name changes are not retained after Compass restarts. Report and export tables show empty well names (21677)
 33. Standards curve graph tab does not appear after an undo-redo action (CSW-72)
Solution: Standards curve tab can be restored by opening analysis setting and clicking on ok button.
 34. In graphs view, when 'stack plots' view is on, consecutive zooming in multi-lane view resets initial zoom action on the Y-axis to display default scale. (CSW-289)
 35. Compass crashes with an internal error dialog when zoom action is performed repeatedly in the graph view in an empty graph area (CSW-313)
Solution: Close the error dialog and launch Compass for SW again, there is no impact on run result data or analysis settings due to this issue.
 36. Export of analysis settings prompts to save in a path different from the run file path (CSW-317)
Solution: One can navigate to intended path and save the analysis settings
 37. Independent Probe Scale option is checked in Lane View even after the setting was unchecked in Graph View (CSW-321)
 38. In images view, capillary selection on Probe 2 images disappears when navigated between standards and samples views (CSW-325)
 39. History log incorrectly displays standard curve name without channel information although peak name assigned to standard curve is not changed (CSW-583)

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

1. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)
2. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272)
3. Cannot copy text to Windows clipboard from Notes or Labels (21289)
4. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
5. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
6. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
7. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
8. Underlines are not aligned for angled and wrapped text Lane labels (21353)
9. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)
10. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)
11. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)

12. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly

Solution: Perform Re-crop to update the Figure Band Label displays (21440)

13. For RePlex runs when “Independent Probe Scale” setting is Off and the user modifies the Contrast in the Lane View, the newly created Figures don’t have the same contrast

Solution: Use Custom Contrast in Crop Regions to adjust contrast as desired (21455)

RePlex™

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
3. Unable to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected.

Solution: Temporarily unselect the Total Protein channel, then the context menu appears, allowing the user to make any necessary changes (19434)

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability

Solution: Use the same fixed contrast levels for added runs (16226)

3. Lane order resets when user switches View from Single to All (21270)
4. Lane order resets after user changes Lane Options (21271)

History

1. Lane View Options and Graph view option changes are not saved correctly in History log (21443)
Solution: Pre-existing access-controlled Jess non-Replex assays that were created with versions earlier than 6.1 will have this issue. If history log entries are critical to the system, perform an initial save of the assay file. Logs of any changes after the first save will be displayed correctly.
2. Removal of Peak name groups entries are not logged (21692).

6.1.0 Compass for SW

Released November 2021

Features

1. Support for the new Stellar assays
2. New image analysis option - High Dynamic Range (HDR) for Jess fluorescence images

Note: The HDR method uses information from multiple exposures to achieve good signal-to-noise for low protein concentration while simultaneously maintaining signal monotonicity at high protein concentration.

3. Jess and Abby RePlex assays are now supported with Access Control Enabled.
4. Saturation warning in experiment tables and red pixel overlay on the Image view when 'Show All Images' is enabled.
5. Support to display both 10-Digit and 11-Digit plate serial number of the assays in the run summary.

Note: When new plates with 11-digit barcode are used with Compass SW 6.0 or older version of software, 'Confirm plate type matches assay' informational message will appear in Jess Start Run dialog window.

Improvements and Fixes

General

1. Updated compass to prevent saving an assay file with invalid well locations (19076)
2. Wes Assay Template Sample Name changes are saved properly (21656)
3. Assay Copy Template - fixed too much white space below the template (21573)
4. Fixed issue - Default Analysis View Options should not be changed for the runs (21447)
Note: Default Analysis View Options are now in sync with assays, not the ones changed for the runs
5. Updated immobilization step for Jess assays to improve instrument performance (21643)
Note: When using existing Jess or Wes assay files in Jess instrument, user will now see a prompt to confirm assay change before performing a run using it (21650).
6. Updated compass to ensure changing exposure for one channel does not impact the lane contrasts of other channels (21548)
7. When Protein Normalization channel is enabled and Normalization is turned Off, PN and PN (%) numbers are shown (21449)
8. In Image view, enabled image "Copy" context option (21360)
9. Change default S/N methods to "Signal-to-Noise Ratio" for Jess default assays (21522).
Note: S/N method "Peak Score (v1-v5)" is renamed to "S/N Compass v5.0" (21659)
10. Enabled turning on/off the password protection of Run Report PDF files, default is on (21579)
11. Abby: Updated to ensure separation movie displays all caps during separation step (21500)
12. Fixed run completed time display issue - run completion time now includes cleanup cycle end time (21538)
13. Fixed the intermittent 'Null Pointer Exception' and 'Graphic is Disposed' internal error dialog issues when playing movie after switching between assay and run views while the instrument is running (21701, 20795)
14. Fixed the issue with Lane view image display - In Lane view of Replex TP runs, when only 'TP Channel' is selected, Lane view image doesn't update to show selected caps when multiple caps are selected (21834)

History

1. Enabled logging of Lane contrast changes in analysis dialog (21108)

2. Apply Settings action is added in History for new Peak Name group for Wes run files (21380)
3. Additional Graph and Lane View options logged in History (21338)
4. Improved image exposure history log entries (21399, 21403, 19056)
5. Log lane contrast setting changes – fixed values and slider levels (19086)

Known Issues

General

1. IV Plot title for Jess run displays "Cycle 1", instead of being blank (18962)
2. History Log - Manual standard changes log message is not easily understandable. (18882)
3. Add Run - Scale incorrect when run files of different sizes are added. (18915)
4. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
5. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
6. Analysis Settings for Peak Names -> Standard Curves -> Concentration - units of precision do not match report. (18905)
7. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
8. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
9. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)
10. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
11. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
12. Start Assay and Blot dialogs can be displayed at the same time. (17467)
13. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
14. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
15. Standard Curve tab disappears upon reanalysis of data if user selects to Cancel settings dialog
Solution: Go to Edit > Analysis... > Standard Curve > OK (19012)
16. Assay Template changes made during run are not saved after the run completes (19050)
17. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
18. Zoom out only works properly for first standard curve graph (21039)
19. Signal to Noise settings – validation checks do not exist to ensure Start < End (21229)
20. File > Close All does not close all run files if one of them needs to be saved (21320)
21. Standard Curve tab disappears after user Adds Run without Std Curve and reanalysis is triggered on second run (21322)
22. Channel Column appears for Peak Groups in Abby run file Reports (21341)
23. Standards Graphs - Exposures in legends should be in black when the plots are stacked (21351)
24. Lane selection for RePlex – Probe 2 lanes not selected if channel added after capillary selection done in Lane View (21388)

25. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (21394)
26. Lane View – Probe 2 Ladder lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (21410)
27. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
28. When Exposure settings contain decimal values, folder name created from 'Export Tables' action in Compass does not display decimals (21613).
29. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)
30. Corrected area column is missing when Normalization is enabled and when ladder cap is set as reference cap (21676)
31. Add runs – 'save', 'close' and 'close all' operations
 - Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed file. The user can then save and close.
 - Selecting File->Close All does not close all the files after saving (19048)
 - When two runs are open, selecting File->Close on a specific run will close the other run(21797).
32. For Locale Chinese and Japanese, in assay perspective well name changes are not retained after compass restarts. Report and export tables show empty well names (21677)

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

1. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)
2. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272)
3. Cannot copy text to Windows clipboard from Notes or Labels (21289)
4. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
5. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
6. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
7. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
8. Underlines are not aligned for angled and wrapped text Lane labels (21353)
9. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)
10. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)

11. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)
12. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly
Solution: Perform Re-crop to update the Figure Band Label displays (21440)
13. For RePlex runs when "Independent Probe Scale" setting is Off and the user modifies the Contrast in the Lane View, the newly created Figures don't have the same contrast
Solution: Use Custom Contrast in Crop Regions to adjust contrast as desired (21455)

RePlex™

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
3. Unable to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected.
Solution: Temporarily unselect the Total Protein channel, then the context menu appears, allowing the user to make any necessary changes (19434)

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability
Solution: Use the same fixed contrast levels for added runs (16226)
3. Lane order resets when user switches View from Single to All (21270)
4. Lane order resets after user changes Lane Options (21271)

History

1. Lane View Options and Graph view option changes are not saved correctly in History log (21443)
Solution: Pre-existing access-controlled Jess non-Replex assays that were created with versions earlier than 6.1 will have this issue. If history log entries are critical to the system, perform an initial save of the assay file. Logs of any changes after the first save will be displayed correctly.
2. Removal of Peak name groups entries are not logged (21692).

2.1 Jess Instrument Software

NOTE: Requires Compass for SW version 6.1 for Stellar assays.

Features

1. Instrument support for Stellar assay (19282)

Improvements and Fixes

1. Self-test improvements – enhanced instrument logs

Known Issues

1. On rare occasions when attempting to close the door, the instrument will not properly recognize the hand sensor input and close (17590)
2. When a run is performed with chemi detection setting other than HDR in the detection profile of Jess assay, chemi channel will not show any data (ESW-142).

Solution: Select High Dynamic Range (HDR) as the Chemi Detection profile to perform chemi based runs on Jess. Details of HDR can be found in Compass for SW user guide.

6.0.0 Compass for SW

Released April 2021

NOTE: If you are using Peggy Sue, Sally Sue or NP1000, Compass for Simple Western version 6.0 and higher does not support these instruments. You can still use software versions 5.x and lower to create, load and run assays on these instruments, and to analyze your data.

Features

1. Support for the new Abby instrument
2. New Lane Annotation tools: Create annotated figures of the Compass Lane view for use in presentations and publications, saving you the hours it would take to manually annotate data using other programs.
3. Analysis perspective **Graph Options** and **Lane Options** selections are now saved with the run file, so that user preferences are saved for the next time the run file is opened (19314)
4. Change default Lane View to not show baseline corrected data. Additionally, move **Baseline On/Off** to Lane Options to support ability to save options with run file (19310, 19566)
5. Default Analysis View options: Graph and Lane View options are saved, created an interface for the default settings, under Edit -> Default Analysis View ... (20742)
6. Support new S/N algorithm improvements, with ability to use existing Compass for SW Peak Score algorithm in versions 5.x and earlier (19289)

Improvements and Fixes

General

1. Improved performance of Lane View display, now with faster rendering of images and overall responsiveness for user interactions
2. Analysis Setting for Peak Names - Loading Controls does not match report (18903)
3. Help menu - Check for Updates, Export Logs and Send Run File options should be disabled during run (18913)
4. Ladder cap should be excluded in Analysis Options -> Peak Names (18944)
5. Capillaries table: Peak Names that are not assigned to Loading control should not be highlighted when Corr. Area is selected (18950)
6. Capillary Groups will not display Rows & Channels for empty groups. (18967, 18969)
7. Fixed display issue of overlapping peak labels for peaks at same MW in different probes (19621)
8. Fixed issue where user could not properly save the run file when using File > Save As... on a file that has the Read-Only attribute set (20806)
9. Standards Graphs - legends are all black for overlay/stacked graphs - should be colored per data trace color (20808)

RePlex™

1. RePlex runs with Total Protein should not display Peak data for Probe 2 Ladder capillary (21195)
2. RePlex runs with Total Protein should not export Primary in Peaks table (21196)
3. RePlex Probe 2 trace colors in report do not match the trace color in Compass (19376)
4. Graph - IR Probe 2 default color is the same as NIR Probe 1 (18993)
5. Peaks table displays Probe 2 Total Protein data for ladder cap (18966)
6. Export: %Total is displayed as 0.0 in Compass for Probe 2 data but in export it is empty (21197)
7. Resolved issue with an incorrect baseline for RePlex Chemi with Total Protein run files. This was an error introduced with the Hot Fix for 19257 in version 5.0.1 (19607)

History

1. History Log does not display old and new values when contrast is changed in Analysis Settings (Edit > Analysis > Lane Contrast) (21108)

Known Issues

General

1. IV Plot title for Jess run displays "Cycle 1", instead of being blank (18962)
2. History Log - Manual standard changes log message is not easily understandable. (18882)
3. Add Run - Scale incorrect when run files of different sizes are added. (18915)
4. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
5. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
6. Analysis Settings for Peak Names -> Standard Curves -> Concentration - units of precision do not match report. (18905)

7. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
8. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
9. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)
10. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
11. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
12. Start Assay and Blot dialogs can be displayed at the same time. (17467)
13. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
14. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
15. When Protein Normalization channel is enabled and Normalization is turned Off, PN and PN (%) numbers are shown (21449)
16. Standard Curve tab disappears upon reanalysis of data if user selects to Cancel settings dialog
Solution: Go to Edit > Analysis... > Standard Curve > OK (19012)
17. Assay Template changes made during run are not saved after the run completes (19050)
18. Should not allow assay Save/Save As with invalid locations (19076)
19. 'Internal Error: Graphic is Disposed' error displayed when attempting to play Separation movie during an Abby run (20795)
20. Plot labels are not displayed correctly (capillary and exposures) when Apply Colors to stacked plots graph options is selected in Preferences for any multichannel runs with RePlex (20835)
21. Zoom out only works properly for first standard curve graph (21039, 21412)
22. Signal to Noise settings – validation checks do not exist to ensure Start < End (21229)
23. File > Close All does not close all run files if one of them needs to be saved (21320)
24. Standard Curve tab disappears after user Adds Run without Std Curve and reanalysis is triggered on second run (21322)
25. Channel Column appears for Peak Groups in Abby run file Reports (21341)
26. Standards Graphs - Exposures in legends should be in black when the plots are stacked (21351)
27. Image View Copy context menu option is disabled for Jess/Abby run files (21360)
28. Lane selection for RePlex – Probe 2 lanes not selected if channel added after capillary selection done in Lane View (21388)
29. Plot label color should be black for stacked plots when "Apply colors to stacked plots" is off (21394)
30. Default Analysis View Options UI display becomes out of sync after View Options are saved in the run file. This is a UI display issue ONLY, as the View options in the Assay saved with the run (*.cbz) are correct (21447)
31. Lane View – Probe 2 Ladder lane location moves when user switches from "View All" to "View Selected"
Solution: User can manually move the ladder lane to desired location (21410)

32. File Open Recent List - file path is not correctly saved for folder name contains comma (21506)
33. German language: Decimal point is not replaced by comma in export table (21673) and in the analysis settings fields of PDF report (21681)

34. Add runs – ‘save’ and ‘close all’ operations

- Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed one. The user can then save and close.
- Selecting File->Close All does not close all the files after saving (19048)

Lane Annotations

NOTE: Software verification only performed on a limited set of font styles and sizes. Some Windows fonts and large sizes will not display well, as with any other free form editing software. Certain resolution size and scaling may also produce undesirable results, use Windows Recommended settings where possible.

1. Custom Band Labels and MW Ladder can overlap when on same side of Figure image (21094)
2. Lane/Band Labels and Notes can overlap for some fonts/sizes (21272)
3. Cannot copy text to Windows clipboard from Notes or Labels (21289)
4. Band Label Linked To field not updated properly after user modifies Peak Name settings (Group name, Peak name, MW) (21299)
5. Lane labels and Caps list incorrect if user hides capillaries in Figure after being created (21316, 21317)
6. Lane Labels are not at the center of the lanes when rotated 45 degree (21327)
7. Default locations for Lane Labels are in reversed order, comparing to Lane view (21348)
8. Underlines are not aligned for angled and wrapped text Lane labels (21353)
9. When adding new MW Region, sometimes the existing Regions will change if the software determines no empty region is available (21390)
10. Exported .tiff image is roughly 50% smaller than .png or .jpg when copied into Word/PowerPoint/etc. (21391)
11. In Notes section, cursor jumps to the beginning of the Notes after "#cbz" button is clicked (21454)
12. When the ladder probe is changed by the user in Analysis Settings, Lane capillary is updated but the Band Labels are not displayed properly
Solution: Perform Re-crop to update the Figure Band Label displays (21440)
13. For RePlex runs when “Independent Probe Scale” setting is Off and the user modifies the Contrast in the Lane View, the newly created Figures don’t have the same contrast
Solution: Use Custom Contrast in Crop Regions to adjust contrast as desired (21455)

RePlex™

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
3. Unable to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected

Solution: Temporarily unselect the Total Protein channel, then the context menu appears, allowing the user to make any necessary changes (19434)

4. In some Re-Plex standards run, Y-axis displays chemi units instead of Fluorescence (21429)

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability

Solution: Use the same fixed contrast levels for added runs (16226)

3. Lane order resets when user switches View from Single to All (21270)
4. Lane order resets after user changes Lane Options (21271)

History

1. Apply Settings not included in History for new Peak Name group for Wes run files (21380)
2. Additional Graph and Lane View options logged in History (21338)
3. In some cases, History does not properly show changes for Image exposure selection for Probe 2 (21399, 21403)

1.3 Abby Instrument Software

NOTE: Requires Compass for SW version 6.0 or higher

Features

1. Support for new Abby instrument

Known Issues

1. None

5.0.1 Compass for SW Released October 6, 2020

Improvements and Fixes

General

1. Fixed issue where calculated Molecular Weight values are incorrect for peaks detected in Probe 1 and Probe 2 in assays where ladder detection is set to Probe 2 (19257)

Known Issues

1. When special characters are used in run file name, Peggy Sue will not play separation movie correctly (21653)

Solution: Do not use special characters in run file name

5.0.0 Compass for SW Released July 21, 2020

Features

1. Support for the new RePlex™ feature on Jess, providing the ability to run two immunoassays within the same capillary to get all your rich protein characterization data from just one sample!
 - Quantify expressed phosphorylated target and total target levels
 - Normalize your data with total protein expression data in the same capillary
 - Greater flexibility in immunoprobe selection
 - More data points per sample
 - Save time and money on consumables

NOTE: RePlex™ assays are not supported with Access Control enabled.

Improvements and Fixes

General

1. Capillaries table displays the aligned position value for Standard Peak 1 position, which is slightly different from the position index value displayed in the Peaks table (17605)
2. Sample fitted peaks for Run Report incorrectly display when user selects Baseline Fit option in View Menu for graphs (16986)
3. Fitted peaks are not displayed in Run Report, even though Fitted Peak option is selected (17734)
4. For Jess assays, automatically expand the Detection node by default to help the user review exposure and ladder settings (18122)
5. Exclude Ladder lane from contrast adjustment to accommodate where user was unable to see ladder peaks in Lane View when sample peaks are 30 times higher or lower (4566)
6. Resolved issue where Graph would show peak values when not selected in Graph Options (9825)
7. Removed Diagnostic option in View Menu for Capillaries table (18509)
8. Fixed issue where same peak label is displayed twice in Graph view (18441)
9. Assay Protocol for Wes Total Protein - "Primary Antibody Time" should be named as "Antibody Diluent Time" (17687)
10. Resolved issue where Standard Curve becomes empty when uses switches from Samples to Standards view (16055)
11. Resolved an issue where the PN Sample peaks corrected area values would change when Protein Normalization is disabled. This was incorrect, the value for these peaks do not change. This will not affect data when using the Protein Normalization feature in Jess (18805)
12. Analysis Settings for Peak Names - Loading Controls will present previously enabled loading controls as enabled in non-selected peak groups. (18938, 18986)

13. For Peggy Sue/Sally Sue/NP1000 assays, display warning message when sample load time is 0 in the assay protocol (18223)
14. Name Peak > New option should exclude ladder capillary (17953)

CFR

1. Audit Trail – improved message for Send Run to include same information as Send Logs (e.g. send_runfile | Uploaded run for [SN], [path] [*.cbz]) (18236)

Known Issues

General

1. IV Plot title for Jess non-RePlex run is "Cycle 1", not blank. (18962)
2. History Log - Manual standard changes log message is not easily understandable. (18882)
3. Add Run - Scale incorrect when run files of different sizes are added. (18915)
4. Report will include legacy "Registration" Standards, although they are not used in analysis, nor are they visible in the Analysis Settings - Standards for Jess or Wes. (18836)
5. If the user creates a name of the form Cn (such as C1) in the assay template, that name cannot be used in the Caps filter for Peak Names in Analysis Options (18753)
6. Analysis Settings for Peak Names -> Standard Curves -> Concentration - units of precision do not match report. (18905)
7. Analysis Settings for Peak Fit - Apply Override range changes are not properly applied to selected caps. (19024)
8. It is possible for Peaks to be improperly labelled if user adds more manual peaks than there are ladder peaks. (18050)
9. It is possible for the check-mark next to a ladder cap to not be removed when ladder peak changes are removed (18996)
10. When trying to force a small/non-existent peak as a ladder peak, analysis will sometimes have issues finding the previous Ladder peaks
Solution: Select Clear from context menu to remove the manual peak (19109)
11. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards. (16351)
12. Start Assay and Blot dialogs can be displayed at the same time. (17467)
13. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files. (16850)
14. ANDI export: In certain cases, peak positions may show slight difference between Compass for SW and Empower/Chromeleon due to peak detect/find algorithms with different software packages used for analysis (18873, 16048)
15. Add runs – 'save' and 'close all' operations
 - Save button will not be enabled for Add Runs when only one file is modified (19042)
Solution: Close the run file that has not been changed, this enables save button for the changed one. The user can then save and close.
 - Selecting File->Close All does not close all the files after saving (19048)

RePlex™

1. Plot colors for RePlex runs may not be in the order expected. (18943)
2. For RePlex assay data, Loading Controls can be used to correct the peaks of the same capillary, for both Probe 1 and Probe 2. When multiple Loading Controls are defined across Probes, corrected area results for some capillaries can be incorrect (18989)
3. Unable to make manual ladder changes for ladder capillary for RePlex assays with Total Protein when all channels are selected

Solution: Temporarily unselect the Total Protein channel, then the context menu appears, allowing the user to make any necessary changes (19434)

Lane

1. Named Peaks are not displayed in Lane View when viewing multiple run files, even if the Named Peaks option in Lane Label dialog is selected. (16227)
2. Contrast enhancement tool in Lane View is currently implemented for single run file only, does not work as expected when viewing data from multiple run files via Add Run capability

Solution: Use the same fixed contrast levels for added runs (16226)

2.0 Jess Instrument Software

NOTE: Requires Compass for SW version 5.0 for RePlex assays. Compass for SW software versions 4.1.0 and lower cannot process Jess instrument run files from 2.0 embedded software

Features

1. Instrument support for RePlex assay (17721)
2. Create explicit cleanup cycle for RePlex assays, which will include an assay node to specify the polymer location (17723)
3. Add parameterized assay nodes and action blocks/units for RePlex Purge and Biotin Labeling to support ChemilA+TP assay type (17766, 17815)

Improvements and Fixes

1. Increase low vacuum self-test summary data precision to 3 digits (17790)
2. When the door is open at power-on for the instrument, the light output from the High-Power LEDs can be less than expected. To mitigate this issue, the embedded software will now turn the HPLED on to 100% briefly when the instrument door closes (17876)
3. Include self-test module exceptions in log files when running tests (18091, 18107)
4. Save specific and unique filenames for each filter alignment test image to aid in troubleshooting (18367)

Known Issues

1. On rare occasions when attempting to close the door, the instrument will not properly recognize the hand sensor input and close (17590)

4.1.0 Compass for SW Released September 25, 2019

Features

1. Online User Guide and Release Notes (16920)
2. Added feature (Help -> Send Run File) that allows customers to send run files directly to ProteinSimple for Technical Support, utilizing a secure cloud storage location (17392)

Improvements and Fixes

General

1. Improved the Jess Start Run dialog for barcode warnings for plate and cartridge (16597)
2. Compass for SW now prevents multiple users or instances of the software from editing the same run file (*.cbz) (17120, 17537)
3. Resolved issue where user could not print Run Reports (16643)
4. Fixed issue where user modified (add/remove) baseline points and changes are not saved properly for Fluorescent channel data in Jess (16475)
5. Resolved issue where the user could not logout from the application when there is an issue with a user account that is currently logged in (17125)
6. Resolved case when Compass for SW software attempts to save an incomplete run file that has not been properly downloaded from the instrument. For this version, Compass will not allow the user to Save the file when it is incomplete (17124)
7. Compass for SW now prevents the user from overwriting an existing assay or run file via the File->Save As... menu operation (17068)
8. Resolved issue where access to Error logs option in instrument properties dialog is not present when connected to non-Jess instruments. To access these logs, click on Logs in the Instrument-> Properties... menu option, then click 'All Logs' button to access list of all log files (16387)
9. Fixed issue where the Start run dialog would not recognize cartridge and plate barcodes if communication to instrument is lost and re-established (17267)
10. Fixed issue where the Start button in the Start Run dialog is disabled if communication to instrument is lost and re-established (16999)
11. Compass for SW will now detect and prompt user when any default assay file is manually deleted or modified (17047)
12. Fixed issue where Compass would display Null Pointer Exception (NPE) error when forcing Standards for the Ladder capillary (16634)
13. Fixed issue where Standards image capillary locations (yellow box) move when IR data channel is selected on Jess (16617)
14. Fixed dialog display title for '*Assay – Detection Profile*' (16697)

15. Improved consistency with 'Assay – Detection Profile' title to display (Chemi), similar to (NIR) and (IR) (17401)
16. Fixed issue where reports cannot be generated for Peggy Sue, Sally Sue, and NP1000 completed run files with embedded software version 4.0 (17543)
17. Fixed Run Report issue where 4PL standard curves are displayed as Linear (17579)
18. Resolved issue when Ladder Capillary is set to 'None' for Peggy Sue or Wes run and changes made to the Peak Fit View region are not applied to the graph (16851)
NOTE – As the Lane View contrast calculation is based on the Analysis View region, Lane View images may look slightly different when compared to images in version 4.0 (17673)

CFR

1. Improved Authorization Server Audit Trail messages to include full file path for sign_off, save_assayfile, save_runfile, open_assayfile, and open_runfile events (17119)
2. Fixed issue where duplicate start run log entries are listed in Run Summary History (17450)
3. Fixed issue where Compass for SW starts an uncontrolled run file when Access Control is On. This happens if communication to instrument is lost and re-established when Start Run dialog is open (17122)
4. Start button for standard capillary cartridge assay in Start Run dialog is now disabled if a blot cartridge is present in the instrument (17438)
5. Compass for SW now records Export Uncontrolled action in History log and Audit Trail (17121)

Known Issues

1. Capillaries table displays the aligned position value for Standard Peak 1 position, which is slightly different from the position index value displayed in the Peaks table (17605)
2. Electropherogram graph scaling issue when viewing multiple run files and one of the run files is not properly aligned to standards (16351)
3. Start Assay and Blot dialogs can be displayed at the same time (17467)
4. Sample fitted peaks for Run Report incorrectly display when user selects Baseline Fit option in View Menu for graphs (16986)
5. Fitted peaks are not displayed in Run Report, even though Fitted Peak option is selected.
Solution: Check option to un-select and select again (17734)
6. Plot Label display is cut off on left side of the plot when Ladder Capillary is set to 'None' and a single capillary is selected. This only applies for Wes and Peggy Sue run files (16850)
7. Named Peaks are not displayed in Lane View when viewing multiple run files, even though Named Peaks option in Lane Label dialog is selected (16227)
8. Contrast enhancement tool in Lane View does not work as expected when viewing capillary data from multiple run files (16226)

1.2 Jess Instrument Software

Improvements and Fixes

1. Added Fluorescence LED calibration routine (17070)
2. Prevent condition where Compass for SW can start a new Assay run. Jess instrument now ensures cartridge barcode image is completed before starting a new run (16581)
3. Resolved issue where some system log files were not being rotated properly (17067, 17034)
4. The Unregulated Vacuum Test can sometimes fail because a single calibration was being used for two vacuum sensors. Modified to calculate the calibration curve for both sensors independently (17045)
5. Improved reliability and robustness of closing the door for Jess instruments (17109)
6. Modified Lamp PWM range for Borofloat solarization effects on Jess (17110)
7. Fixed issue where calibration could not be performed on instrument the day before the Light Saber tool expires (16633)
8. Added capability to detect background artifacts from cartridge (17269)
9. Display Instrument name in service console browser tab page title for main page (17436)
10. Add log entry in embedded.log file for r-squared when determining pump curve in instrument high vacuum calibration routine (17434)
11. Modified high vacuum calibration routine to include additional venting and settling time when determining setpoint (17452)
12. Added log entries for any exceptions encountered in Camera Focus Calibration setup (17484)

Known Issues

1. On rare occasions when attempting to close the door, the instrument will not properly recognize the hand sensor input and close (17590)
2. On rare occasions, the Jess instrument will not boot properly after a power cycle, displaying purple on the instrument status light. Contact Technical Support for further troubleshooting and resolution (17040)

4.1 Peggy Sue/Sally Sue/NP1000 Instrument Software

Improvements and Fixes

1. Resolved issue where some system log files were not being rotated properly (17067, 17034)
2. Fixed issue where dark background images were not being cropped (17066, 17492)

4.0.1 Compass for SW Released August 9, 2019

Improvements and Fixes

1. New user guide has correct reference to EMC standard (17614).

4.0.0 Compass for SW Released June 8, 2018

Features

1. Support for Jess instrument – application support for two-color fluorescence detection, protein normalization, expanded self-test, cartridge information tracking, and traditional western blot imaging.
2. Analysis Options panel - Allows for easier access to channel exposure selection and peak naming from the Analysis view without having to open the preferences dialog.
3. Ability to name peaks from Graph/Lane view and Peaks table
4. Improved layout and organization of Analysis/Preferences dialog.
5. Support for multiple Standard Curves.
6. Ability to configure Auto Lock Inactivity timeout in Preferences. Default value is 20 minutes.
7. Display of peak name labels in Lane View
8. New Export logs feature, allowing the user to gather all logs from the instrument and Compass into a single file.
9. Improved grouping and configurability for generating Run Report.
10. Support for encryption of PDF file for Run Reports
11. For advanced users, ability to view multiple channel images.

Improvements and Fixes

1. Improvements to High Dynamic Range analysis - HDR 4.0 handles a wider range of burnout cases. HDR 4.0 is less sensitive to the level at which burnout occurs. When burnout occurred at a low level the classical HDR method had trouble picking the optimal exposure and interpolating between exposures (13882, 14941, 14856, 14963, 9905, 9794)
2. Improved Standards peak find – additional detection and correction of outliers (9469)
3. Peggy Sue/Sally Sue – support Ladder settings per cycle, so that user can have different ladder cap settings for different cycles. Allows user to fix data if the ladders for one or more cycles are bad or missing (4770, 4771)
4. Assay Protocol – removed user ability to modify wash related steps for Wes/Jess assays. This was done to prevent instrument failure when too many washes were specified by the user (10143)
5. Increase pl precision on graph and header to match Peak Table (13884)
6. Adjust Lane View contrast scaling to the current selected view region (8093)
7. Resolved NPE error where user was unable to analyze and display data for a Peggy Sue run with Capillary boxes empty error (10327)
8. Corrected immobilization wash load time for 13 cap assay (13943)
9. Resolved issue where grouping order shown in Peak Groups was not consistent with Cap Groups and Capillary Groups (9975)
10. Resolved issue where Charge Assay - Separation Type Voltage is not supported on Peggy Sue. User would see a Java NPE error upon changing the exposure time (14985)
11. Resolved issue where peak labels and peak results disappear for the ladder cap after adding a ladder peak on graph (14005)

12. Hide "Cycle" column in Experiment pane for single-cycle runs (3700, 4017, 4297, 15100)
13. Added support for fractional time zones (15902)
14. Resolved issue where Compass does not properly display the Instrument date and time after using 'Set to PC Time' function (15965, 15922)

NOTE: 'Check for Updates' feature in Compass software will not update to version 4.0.0. User will need to download installer from ProteinSimple website (<http://proteinsimple.com/compass/downloads/>) and perform a full installation.

3.1.8 Compass for SW Released June 8, 2018

Improvements and Fixes

1. Access Control – resolved issue where login button is enabled after user enters in a value in the password field and then removes the entry (15678)
2. Resolved issue where Lane View label font is corrupted after applying Windows 10, April 2018 update, version 1803 (15690)

3.1.7 Compass for SW Released December 15, 2016

Features

1. New High Dynamic Range (HDR) Detection Profile for size and charge assay (Wes, Peggy Sue, Sally Sue)

The Detection profile is set as default and can be disabled from the Assay protocol and replaces the old Multi-Image Analysis. For Wes, the higher dynamic range is achieved by using a more effective algorithm to create the High Dynamic Range multi-image analysis, refresh substrate from the assay plate prior to each exposure and minimizing the time between substrate aspiration and Image. No substrate refresh takes place in Peggy Sue and Sally Sue.

For Wes, the embedded or instrument software must be updated to v2.5 or higher.

Old runs (performed with Compass for Simple Western v3.0 or Compass v2.7 or older) can be analyzed with the new algorithm for High Dynamic Range.

The substrate refresh requires that 15 μ L of substrate instead of 10 μ L be loaded /well for all Wes assays. A window with a warning to check for correct volume pops up after the Start run button is pressed.

3.0 Compass for SW Released August 15, 2016

Features

1. New name and logo

Old = Compass 

New = Compass for Simple Western 

2. Support for Low MW assay

- a. New assays for Low MW 2-40 kDa are added for Wes, Peggy Sue, and Sally Sue.
- b. Split buffer choice eliminated since it's the default buffer for all assays.

NOTE: Embedded update to v3.4.29684 required for Peggy Sue and Sally Sue to run new Low MW assays

NOTE: Low MW runs generated by v3.0 cannot be opened with Compass v2.7.

3. Multiple Exposure View

Turn on "All Exposures" view from **Graph Options** - checking this option will overlay all the chemiluminescence exposure times acquired for a particular run. The peak that corresponds to the exposure time selected in the Analysis window under Images will be outlined in blue and the Peaks table below the graph window displays the data corresponding to the selected exposure time

4. Run Reports - run report in PDF

- a. New menu item – File > Run Report...
- b. Run Report dialog – options for report sections and report name
- c. A PDF report is created with the options specified.

5. History Details - Old and new values changed in Analysis Settings are logged

6. Command Log - Added "Save File As..." button to save the log file

7. Peak Fit View Options

- a. Added a "View" setting to control the default view to Full or Analysis.
- b. Added an "Apply" button to see analysis changes without closing the dialog.
- c. Removed the "Restore Original" button.

8. Grouping Options Moved to Preferences - Simplified the Grouping command to toggle the Grouping display without a dialog

9. Lane Contrast in Analysis window - Edit>Analysis>Lane Contrast

Check the radio button for Fixed Level. The numbers displayed refer to the current White and Black levels for the run being analyzed. Change these settings and click Apply to see how they impact the contrast on the lane view. Once satisfied with the changes, click OK. When the you select Fixed Level, the slider tool is no longer available in the lane view. Lane Contrast settings from the assay or run analysis windows can be exported to be applied to other assays or runs.

NOTE: Files created or saved with the Lane Contrast feature locked cannot be opened by v2.7.

10. Raw Baseline

Improvement in how the raw baseline is calculated. Raw baseline for the shortest exposures is shifted (typically lower than previously observed if you are looking at a data obtained with a prior version). That said, the peak area for all exposures will remain the same and will not be affected by the shift of the raw baseline. Area to baseline ratio, on the other hand will change from previously calculated with earlier versions of the software.

11. Other Misc. Changes

- a. Allow size analysis without ladder peaks. Standards for fitting will be used for MW.
- b. Cleaned up peak table so only Standards or Registration peaks are displayed in peak table.
- c. Reduced the plot line width of exported graph.
- d. Renamed the view menu items and changed the corresponding icons on the toolbar.
 - i. View > Multiple View has been renamed as View > All
 - ii. View > Single View has been renamed as View > Selected
- e. Added a Self Test Confirmation dialog.
- f. Added a "Save File As" button in Command Log dialog to save the log.
- g. Added the instrument software version to Comment in Command Log after software upgrade

2.7 Compass Software Released August 31, 2015

Features

1. Support for Split Running Buffer assays for Peggy Sue and Sally Sue and Wes size assays. Peggy Sue and Sally Sue also require embedded software v3.3 or newer
 - a. "Split Running Buffer" will be displayed on the column header for split running buffer assays.
 - b. "Upper RB" and "Lower RB" will be displayed in the Start Wizard for split running buffer assays.
 - c. "Split Running Buffer" will be displayed in Kit Info for the runs with split running buffer assays.
 - d. "Split Running Buffer" will be appended to the Plate S/N Info for Wes runs
2. Export and import assay template now supports *.csv and *.txt formats
3. Support copy and paste assay well contents from Excel and between Compass windows, including both names and attributes
4. Fixed bug to make it possible to create a view for Group Plot if the group statistics for one of the groups is zero
5. Fixed bug that created Duplicate Peak Name columns in Capillaries table when there are multiple identical peak names

2.6.8 Compass Software *Released July 24, 2015*

Features

1. Wes plate barcode is captured for runs with embedded software v2.2.25328 and newer. Plate S/N number is displayed after run is complete in the Run Status tab.
2. View Region – Lower and Upper values now reflect the option selected (Analysis, Full, or Custom). Previously only custom values are displayed.
3. Download and open last completed run if Compass was disconnected during the run.
4. Fixed bug that caused manual standards to be lost after reanalysis, if Analysis dialog was cancelled previously.

2.6.7 Compass Software *Released May 31, 2015*

Features

1. Support Pipet Wash for Sally Sue and Peggy Sue instruments, with instrument software v3.2 and up.
2. This release changes the default Size and Total Protein assays for Sally Sue and Peggy Sue to be compatible with Running Buffer 2 Rev B (Size and Total Protein assays 12-230 kDa). When File>New is selected, Rev. B assays will be created.

2.6 Compass Software *Released September 15, 2014*

Features

1. Support for new Total Protein Size assays.
2. Support for new 66-440 kDa Size assays.
3. Support for updating instrument software from Compass

2.5 Compass Software *Released January 31, 2014*

Features

1. Support for new instrument, Wes
2. Support for new instrument, Sally Sue
3. Support for new instrument, Peggy Sue
4. New Loading and System Control functionality
5. New standard curve fit functionality

2.4 Compass Software *Released June 28, 2013*

1. Support for 21 CFR Part 11

2.3 Compass Software *Released November 30, 2012*

1. Support for new instrument, Peggy

2.2 Compass Software *Released June 30, 2012*

1. Support for new instrument, Sally
2. New grouping and statistics functionality

2.1 Compass Software *Released November 30, 2011*

1. Lane view improvement for sizing data

2.0 Compass Software *Released October 31, 2011*

1. Support for new instrument, Simon.